


# MICROVUE™ Creatinine Assay Kit

Bone Health

An enzyme immunoassay for the  
quantitation of Creatinine in urine

For Research Use Only.  
Not for use in diagnostic procedures.

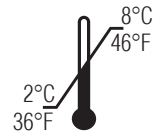

**QUIDEL**®  
CORPORATION  
SPECIALTY PRODUCTS  
RESEARCH TO RAPIDS®

REF 8009

# MicroVue™ Creatinine Assay Summary

## Reagents and Samples Preparation

- Prepare Color Solution (*add 1 mL Stop Solution to each required vial*)
- Prepare urine samples according to PI instructions and dilute 1:40 with deionized or distilled water (*e.g. 20  $\mu$ L sample + 780  $\mu$ L water*)
- Dilute Standards and Controls 1:40 with deionized or distilled water (*e.g. 20  $\mu$ L sample + 780  $\mu$ L water*)

## Assay Procedure

Pipette **50  $\mu$ L** of each diluted Standard, Control, or urine into assay wells (*Use deionized water as the "0" standard*)



Carefully pipette **150  $\mu$ L** of Working Color Solution into assay wells



Incubate **30  $\pm$  2.5 min** at 18 – 28°C



Read the Optical Density at 490 nm  
Analyze the assay results using a linear curve fit  
( $y = mx + b$ )

## SUMMARY AND EXPLANATION

Creatinine is a molecule with a molecular weight of 113.1 g/mol. Creatinine is found in muscle tissues and is excreted into the circulation at a constant rate. It is removed from the plasma by glomerular filtration in the kidney and excreted into the urine.

## PRINCIPLE OF THE PROCEDURE

The MicroVue Creatinine assay is a quantitative, colorimetric assay based on a modified Jaffe method where alkaline picrate forms a colored solution in the presence of creatinine.

## REAGENTS AND MATERIALS PROVIDED

### 96 Assays for the Creatinine assay

The MicroVue Creatinine assay contains the following:

<b>A Creatinine Standards</b>	<b>Parts 4225 – 4227</b>	<b>0.3 mL each</b>
<b>B A = 5 mmol/L, B = 20 mmol/L, C = 40 mmol/L</b>		
<b>C Liquid Creatinine in phosphoric acid</b>		
<b>L Low Controls</b>	<b>Parts 4228</b>	<b>0.3 mL each</b>
Liquid Creatinine in phosphoric acid		
<b>H High Controls</b>	<b>Parts 4229</b>	<b>0.3 mL each</b>
Liquid Creatinine in phosphoric acid		
<b>① Color Reagent</b>	<b>Part 4224</b>	<b>3 x 7 mL</b>
Picric acid (0.14%) in sodium borate and SDS		
<b>② Stop Solution</b>	<b>Part 4960</b>	<b>15 mL</b>
1N NaOH		
<b>Microassay Plate</b>	<b>Part 0673</b>	<b>12 each</b>
Eight-well non-coated strips in a frame		

## MATERIALS REQUIRED BUT NOT PROVIDED

- Timer (60 minute range)
- Tubes for dilution of samples, standards, and controls
- Calculator or other computational method to validate the assay
- Clean, unused microassay plates and/or test tubes and racks
- Container for wash buffer dilution
- Wash bottle or other immunoassay washing system
- Adjustable multichannel pipette (8 or 12 channels) or repeating micro-pipettes (optional)
- Clean pipettes, 1 mL, 5 mL, and 10 mL
- Micropipettes and pipette tips
- Plate reader capable of optical density readings between 0.0 and 2.0
- Deionized or distilled water

## WARNINGS AND PRECAUTIONS

1. For Research Use Only. Not for use in diagnostic procedures.
2. Treat specimen samples as potentially biohazardous material. Follow Universal Precautions when handling contents of this kit and any patient samples.
3. Dispose of containers and unused contents in accordance with Federal, State and Local regulatory requirements.
4. Use the supplied reagents as an integral unit prior to the expiration date indicated on the package label.

5. Wear suitable protective clothing, gloves, and eye/face protection when handling contents of this kit.
6. Store assay reagents as indicated.
7. Do not use Strips if pouch is punctured.
8. Test each sample in duplicate.
9. 1N NaOH may be harmful upon ingestion and can cause irritation to the skin or eyes. Avoid contact with skin, eyes or clothing. If contact is made, wash with water. If ingested, call a physician.
10. The MicroVue Creatinine Color Reagent (picric acid) may be harmful upon injection. Avoid contact with skin, eyes, or clothing. If contact is made, wash with water. If ingested, call a physician. Do not let picric acid dry as the crystals are explosive. If this solution is spilled, flush with water and dispose of the material in an appropriate manner.
11. Use of multichannel pipettes or repeat pipettors is recommended to ensure timely delivery of reagents.
12. For accurate measurement of samples, the addition of samples and standards must be precise. Pipette carefully using only calibrated equipment.
13. Perform a standard curve with each assay.
14. Creatinine values from nonhuman species have not been evaluated with the MicroVue Creatinine Kit.
15. This assay is validated for urine samples only.

## **REAGENT PREPARATION**

**Bring all reagents and materials to 18–28°C before use.**

After removing the needed reagents and materials, return the unused items to their appropriate storage temperatures (see *STORAGE*).

Refer to Table 1 for the amounts of reagents and materials required per number of wells.

### **Working Color Solution**

Prepare Working Color Solution by adding 1 mL of Stop Solution to each required bottle of Color Reagent (see Table in *ASSAY PROCEDURE* section). Color Reagent should be at room temperature prior to use. If a precipitate develops, warm the bottle at low temperature (37°C) to dissolve the precipitate, but equilibrate to room temperature before use.

## **STORAGE**

**Store the unopened kit at 2–8°C.**

**All reagents must be brought to 18–28°C before use.**

## **SPECIMEN COLLECTION AND STORAGE**

**Handle and dispose of all specimens using Universal Precautions.**

Keep the urine samples at 2–8°C if testing within 2–3 days of collection. Store at -20°C for longer storage. Do not subject samples to more than four (4) freeze/thaw cycles.

## ASSAY PROCEDURE

Read entire product insert before beginning the assay.

See **WARNINGS AND PRECAUTIONS** and **REAGENT PREPARATION**.

Determine the amount of Color Reagent required for the number of strips to be used.

# of Strips	2	4	6	8	10	12
# of Samples (tested in duplicate)	2	10	18	26	34	42
Color Reagent (bottles)	1	1	1	2*	2*	2*

\*When more than one bottle is used, combine the contents, and mix prior to use.

### Sample Preparation

1. Mix urine samples and allow to settle for 15–20 minutes, or centrifuge at low speed (200 x g) for at least two (2) minutes. This prevents debris from interfering during pipetting.
2. Dilute Standards, Controls, and urine samples 1:40 with deionized or distilled water (e.g. 20  $\mu$ L sample + 780  $\mu$ L water). Use deionized or distilled water as the zero “0” standard.
3. Place desired number of 1 x 8 strips in the Stripwell Frame.
4. Prepare Working Color Solution by adding 1 mL of Stop Solution to each required bottle of Color Reagent (see Table above). Color Reagent should be at room temperature prior to use. If a precipitate develops, warm the bottle at low temperature (37°C) to dissolve the precipitate, but equilibrate to 18–28°C before use.

### Incubation

5. Add 50  $\mu$ L of diluted Standard (0 (water), A, B, C), Control or urine sample to each well of the 1 x 8 strips in the frame.
6. Carefully add 150  $\mu$ L of the Working Color Solution to each well.
7. Incubate  $30 \pm 2.5$  minutes at 18–28°C.

### Read

8. Read the optical density at 490 nm. Assure that no large bubbles are present in wells and that the bottoms of the strips are clean. Read strips within **10 minutes** of completion of the incubation step.
9. Use quantitation software with a linear regression equation to analyze the MicroVue Creatinine assay results.
10. Determine concentration of samples and Controls from the standard curve.
11. Control values should be within the range specified in the Certificate of Analysis supplied with the kit.

### QUALITY CONTROL


The Certificate of Analysis included in this kit is lot specific and is to be used to verify that the results obtained by your laboratory are similar to those obtained at Quidel Corporation. The optical density values are provided and are to be used as a guideline only. The results obtained by your laboratory may differ.

Quality control ranges are provided. The control values are intended to verify the validity of the curve and sample results. Each laboratory should establish its own parameters for acceptable assay limits. If the control values are NOT within your laboratory's acceptance limits, the assay results should be considered questionable and the samples should be repeated.

## INTERPRETATION OF RESULTS

### Representative Standard Curve

Standard Creatinine levels: 0, 5, 20, 40 mmol/L

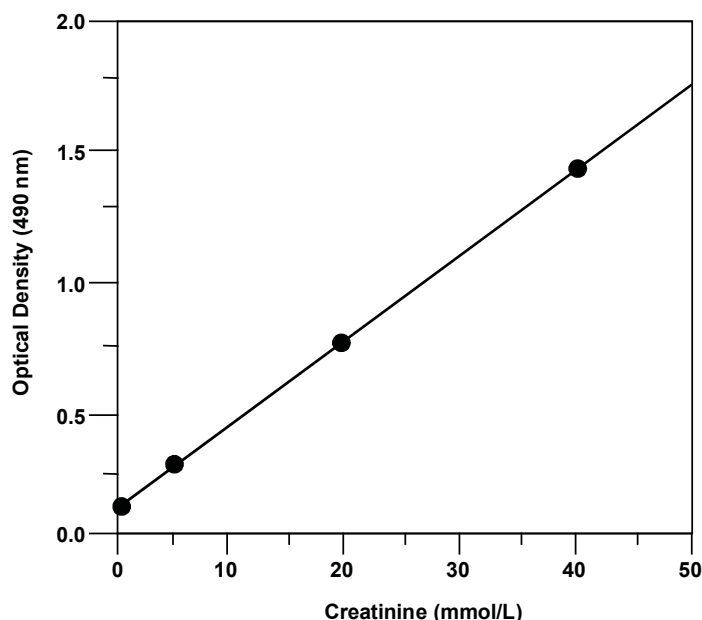

## PERFORMANCE OF THE TEST

### Precision

Within-run and between-run precision were determined by assaying six urine samples in ten different runs. Typical results are provided below.

Creatinine (mmol/L)	Within-run <sup>1</sup> CV (%)	Between-run <sup>2</sup> CV (%)
4.8	1.6	6.9
8.3	2.1	4.6
13.1	1.7	1.6
17.1	1.4	5.4
22.8	2.0	2.0
28.3	1.1	5.4

<sup>1</sup> n = 40 replicates

<sup>2</sup> n = 10 runs

### Recovery – Spike Recovery

Sample	Endogenous (mmol/L)	Added (mmol/L)	Observed (mmol/L)	Recovery (%)
1	3.9	8.4	12.8	106
2	11.0	8.4	20.4	112
3	14.7	8.4	23.8	108


### Recovery – Linearity

Linearity was determined by serially diluting samples and comparing observed values with expected values. Typical results are provided below.

Sample	Dilution Factor	Observed (mmol/L)	Expected (mmol/L)	Recovery (%)
1	neat	36.2	–	–
	1:2	18.4	18.1	102
	1:4	9.1	9.1	100
	1:8	4.5	4.5	100
2	neat	32.2	–	–
	1:2	16.3	16.1	101
	1:4	8.3	8.1	102
	1:8	4.1	4.0	103
3	neat	22.3	–	–
	1:2	11.2	11.2	100
	1:4	5.4	5.6	96
	1:8	2.5	2.8	89

### Accuracy

Accuracy was determined by comparing creatinine values with those assayed by autoanalyzer at a reference laboratory.

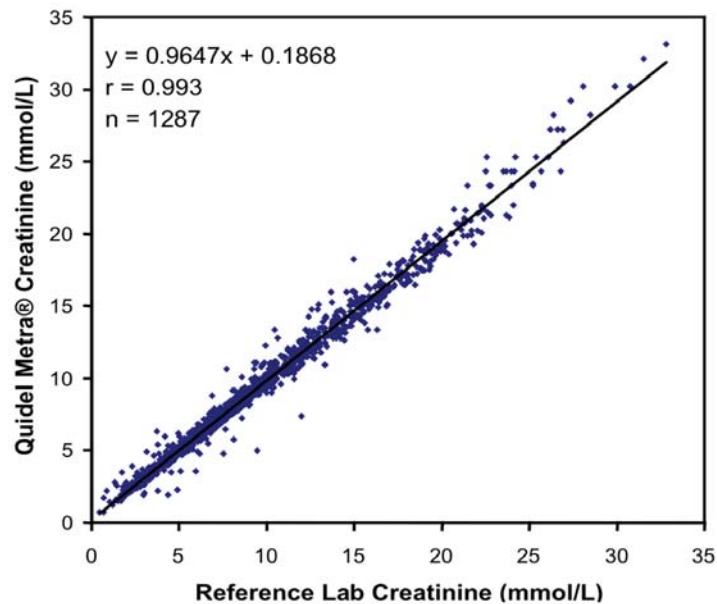

## ASSISTANCE

To place an order or for technical assistance, please contact a Quidel Representative at 800-524-6318 or 408-616-4301, Monday through Friday, between 8:00 a.m. and 5:00 p.m., Pacific Time. Orders may also be placed by fax at 408-616-4310.

For services outside the U.S., please contact your local distributor. Additional information about Quidel and Quidel's products and distributors can be found on our website at [www.quidel.com](http://www.quidel.com).

## REFERENCES

1. Cook JGH. Factors influencing the assay of creatinine. *Ann. Clin. Biochem.* 1975; 12:219-232.
2. Heinegård D, Tiderström G. Determination of serum creatinine by a direct colorimetric method. *Clin. Chim. Acta.* 1973; 43:305-310.
3. Jaffe M. Ueber den neiderschlag, welchen picrinsäure in normalem harn erzeugt und über eine neue reaction des kreatinins. *Z. Physiol. Chem.* 1886; 10:391-400.
4. Owen JA, Iggo B, Scandrett FJ, Stewart CP. The determination of creatinine in plasma or serum, and in urine; a critical examination. *Biochem. J.* 1954; 58:426-437.
5. Tobias GJ, McLaughlin RF, Hopper J. Endogenous creatinine clearance: a valuable clinical test of glomerular filtration and a prognostic guide in chronic renal disease. *N. Engl. J. Med.* 1962; 266:317-323.
6. Centers for Disease Control. Recommendations for prevention of HIV transmission in health-care settings. *MMWR* 1987;36 (suppl no. 2S):001.





Manufacturer



Catalog Number



Consult Instructions for Use



Temperature Limitation



Contents / Contains



Instructions for Use on CDROM



Contains sufficient for <n> tests

---

**REF** 8009 – MicroVue™ Creatinine Assay Kit



Quidel Corporation | 10165 McKellar Court  
San Diego, CA 92121 USA | [www.quidel.com](http://www.quidel.com)

**0279G (2009/03)**