

Technical Information

Anti C4d, FITC

Cat. No.: BI-RC4D-FITC
Volume: 100 µl (125 tests)
Method: **Flow Cytometry**
Sample type: serum
Application:

The antibody has been tested for the flow cytometric detection of cell- or solidphasebound C4 and C4d split product of alloantibody-triggered C4d deposition to FlowPRA® Class I and II Screening Test beads (One Lambda, Inc., Canoga Park, CA, USA). This assay ([C4d]FlowPRA) is applicable for the detection of complement-fixing anti-HLA alloantibodies. For [C4d]FlowPRA (see working protocol). A final dilution of 1:50 is recommended.

Species: Human

Intended use:

C4d is the degradation product of the activated complement factor C4, a component of the classical complement cascade, which is typically initiated by binding of antibodies to specific target molecules. Detection of C4d is regarded as an indirect sign, a “footprint” of an antibody response against the allograft. The majority of publications describe C4d as an important biomarker in kidney transplantation but also in heart, liver, and other transplants.

Intended applications:

- Humoral transplant rejection

For further information please contact / Für weitere Informationen wenden Sie sich bitte an /
Pour plus d'informations, veuillez contacter:

www.tecomedical.com



A EUROBIO SCIENTIFIC COMPANY

Switzerland / Headquarters

TECOmedical AG
Gewerbstrasse 10
4450 Sissach

Phone +41 61 985 81 00
Fax +41 61 985 81 09
Mail info@tecomedical.com

Germany

TECOmedical GmbH
Wasserbreite 57
32257 Bünde

Phone +49 52 23 985 99 99
Fax +49 52 23 985 99 98
Mail info@tecomedical.com

Austria

TECOmedical AG

Phone 0800 20 40 66
Fax 0800 20 40 55
Mail info@tecomedical.com