



# Quick guide to updating the F200 software

## USB drive preparation:

To install the update on the F200, all you need is an empty USB stick that **must be formatted in FAT32** and the corresponding update file.

To format the USB stick in **FAT32**, please connect the USB stick to the PC and then right-click on the USB drive and select "Format"

In the new window, make sure that '**FAT32**' is selected, click on "Start" to format the USB stick to the desired format.

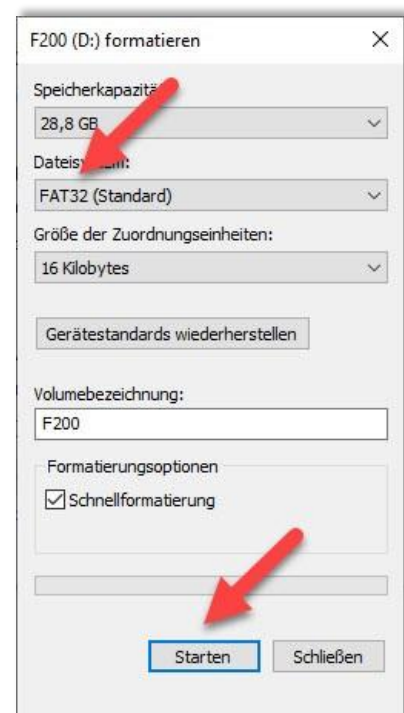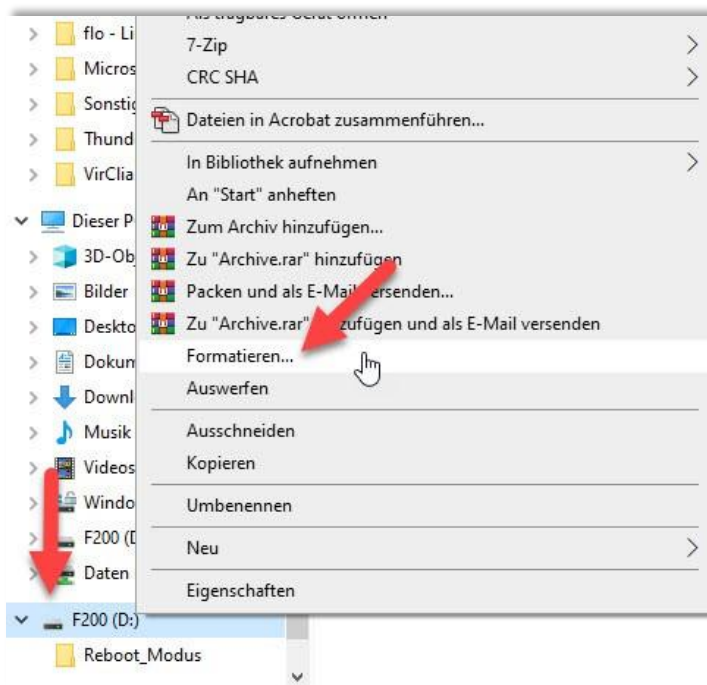


Finally, you only need to copy the file "FluroTest\_V001.067.004.bin" to your USB stick. Your USB stick is now ready for the update.


Procedure on the F200 device:



The prepared USB stick is now inserted into one of the four free slots on the F200 and the device is started.

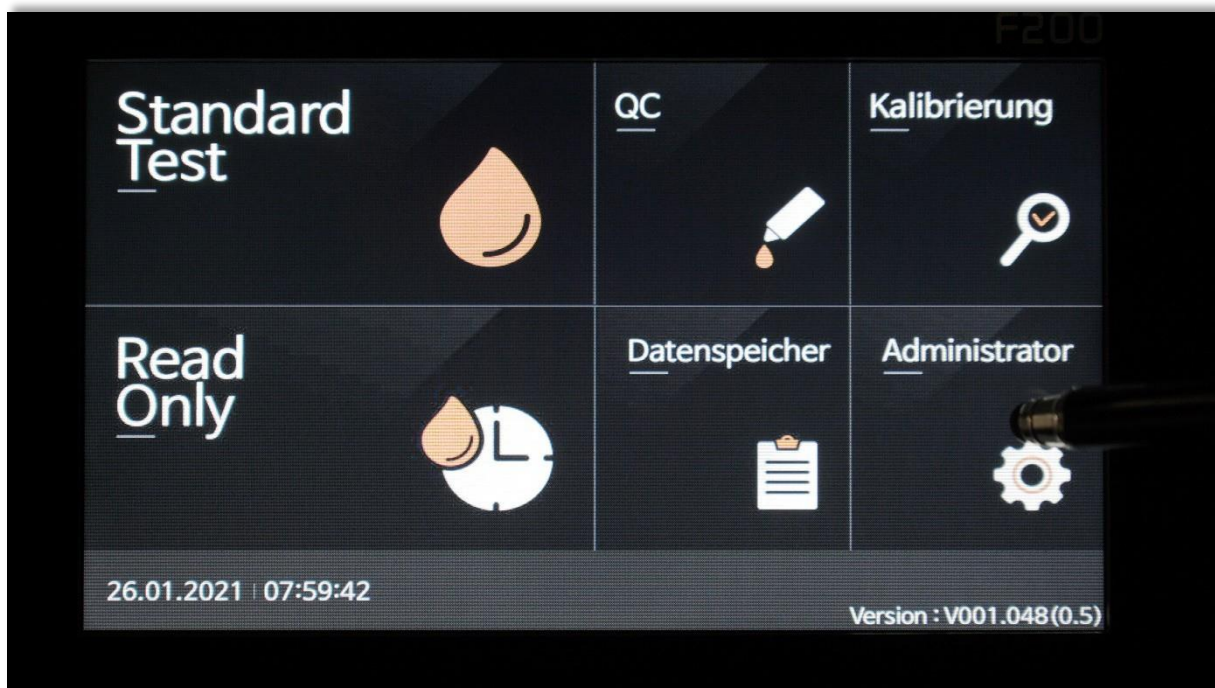

Choose „Administrator“.

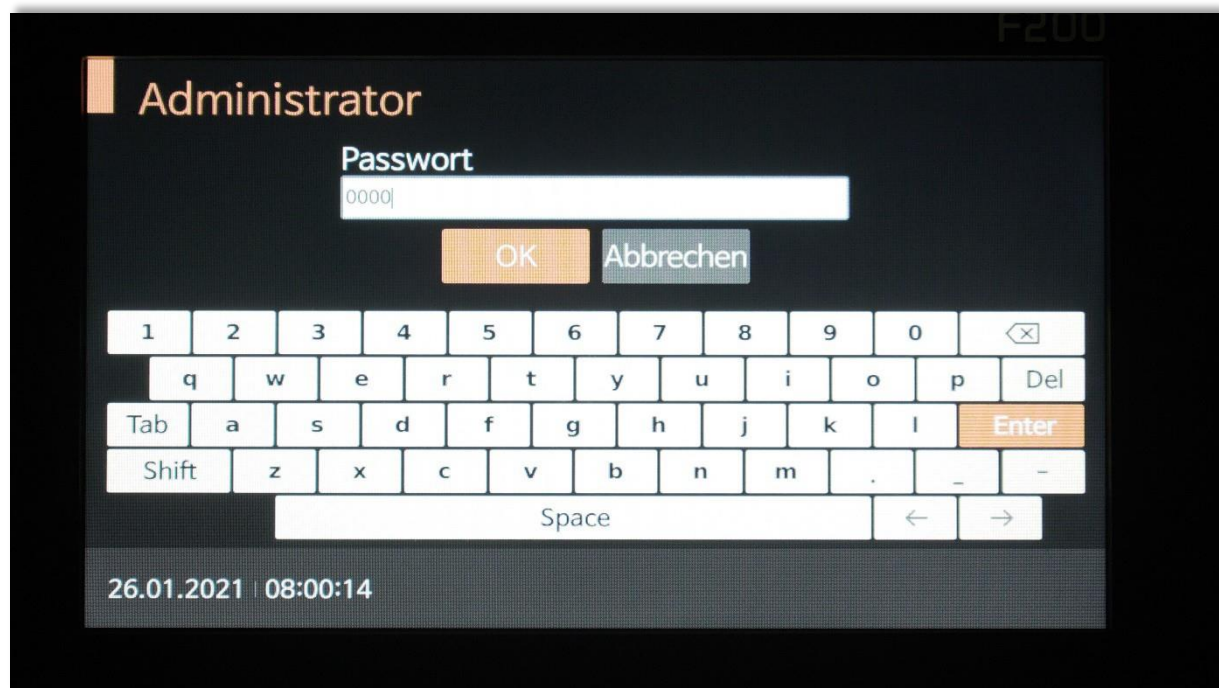


Password: „0000“. Click „OK“.



Click on „update“.



Confirm by clicking „ok“.

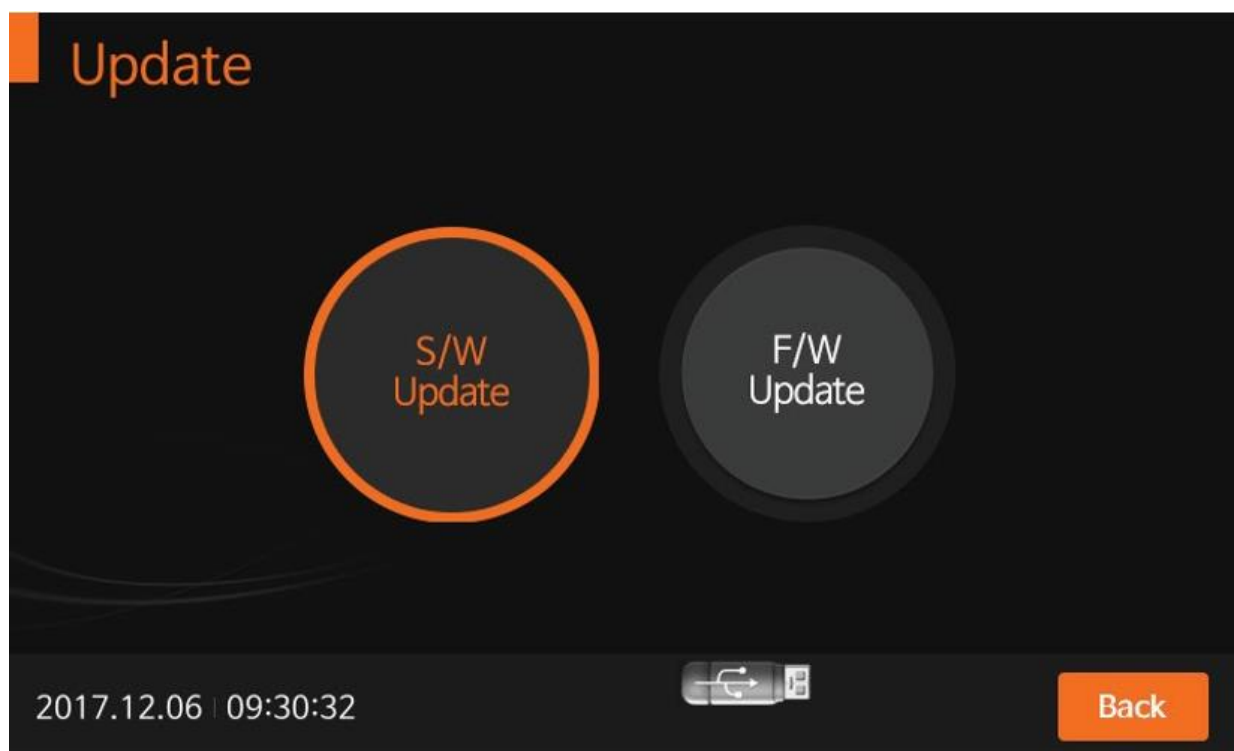

The update is installed after approx. 90 seconds. Please restart the device and remove the USB stick. Your device is now up to date. After the restart, the latest software version is displayed in the bottom right corner of the screen.